
HEAVY-DUTY OBJECT HOLDER

This thing is extremely heavy duty and its primary purpose here is to fixture objects on the HOLOGRAPHIC STEREOGRAM PRINTER when it is used in its object rendering mode, (See the Handout,

Vertical Height Mounting Assembly

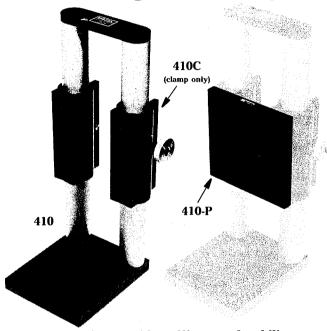


• Dual-rod design provides stiffness and stability

The Model 410 Mounting Assembly positions heavy optical components and assemblies at various elevations with excellent stability. The assembly consists of two 1.5 in. (38 mm) diameter solid steel rods attached to a heavy-duty base and connected together at the top, with an extra-long rod-clamp, Model 410C, on each rod. These clamps have three line-contact pads over the full length of their bore to ensure stable mounting of components. Components can be attached directly to the clamps, or for more versatile mounting of components, the optional Model 410-P Mounting Plate features a standard pattern of tapped holes for-mounting virtually any component.

Vertical Height Mounting Assembly

	English Model	Metric Model
Vertical Height Mounting		
Assembly	410	M-410
Clamp Only	410C	M-410C
Mounting Plate	410-P	M-410-P

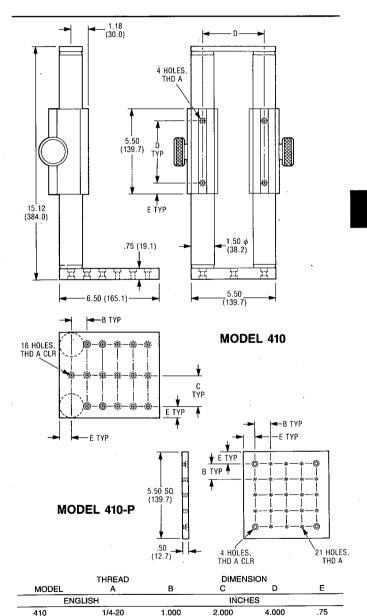

410-P

M-410

M-410-P

1/4-20

METRIC

100.0

.75

199


2.000

MILLIMETERS

1.000


CONVERTING THE HOLOGRAPHIC STEREOGRAM PRINTER TO MAKE HOLOGRAMS OF DIFFUSELY REFLECTING OBJECTS) or on the EXPERIMENTAL TABLE to do the same thing in the SINGLE BEAM TRANSMISSION MASTER HOLOGRAM WITH MIRROR set up. The Model 410-P Mounting Plate is attached to the Model 410C Clamps, and an Enco Model 418-4525 Ground Open End Slotted Angle Plate provides a convenient base for the object to be holographed. Although heavy and stable, (Guess its weight!* Answer below.) it is a good idea to attach the Model 410 to the Tabletop with screws or Newport Model BC-2 Base Clamps.

TO USE, attach the Object to the Holder and loosen the knobs on the Model 410-C Clamps until the Object is at the right height from the vantage point of the Holographic Plateholder. Move the whole unit to get the Object centered horizontally and positioned longitudinally.

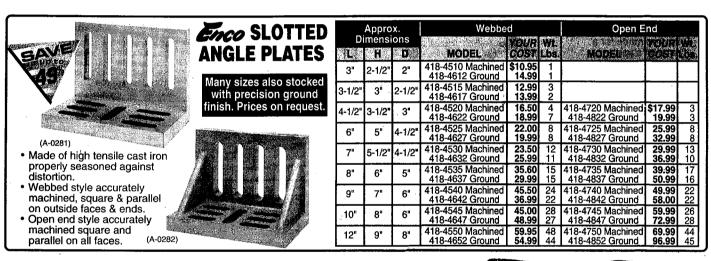

The Slotted Angle Plate can be attached to the Model 410 with 1/4"-20 screws and washers, and oriented up, down or sideways, with Objects hot glued directly to it, or the Objects glued to a Magnetic Base so that they can be easily removed and positioned during a holographic multi-channel exposure.

The latest incarnation of the beast, **Model 410-RC** uses a Siamese twinned set of clamps.

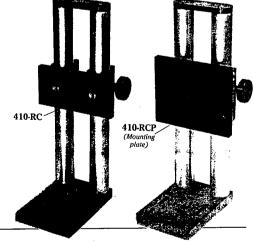

^{*.} According to the scale in the Nurse's Office, the fully-decked out Model 410 plus Right Angle Platform is 27 lbs.